

Solvent Data Chart

	1H Chemical Shift (ppm from TMS) (multiplicity)	(zH) UHľ	¹³ C Chemical Shift (ppm from TMS) (multiplicity)	JCD (Hz)	1H Chemical Shift of HOD (ppm from TMS)	Density at 20°C	Melting point (°C)	Boiling point (°C)	Dielectric Constant	Molecular Weight
Acetic Acid-d ₄	11.65 (1) 2.04 (5)	2.2	178.99 (1) 20.0 (7)	20	11.5	1.12	16.7	118	6.1	64.08
Acetone-d ₆	2.05 (5)	2.2	206.68(1) 29.92 (7)	0.9 19.4	2.8*	0.87	-94	56.5	20.7	64.12
Acetonitrile-d ₃	1.94 (5)	2.5	118.69 (1) 1.39(7)	21	2.1*	0.84	-45	81.6	37.5	44.07
Benzene-d ₆	7.16 (1)		128.39(3)	24.3	0.4	0.95	5.5	80.1	2.3	84.15
Chloroform-d	7.24 (1)		77.23	32.0	1.5*	1.50	-63.5	61-62	4.8	120.38
Cyclohexane-d ₁₂	1.38 (1)		26.43 (5)	19	0.8	0.89	6.47	80.7	2.0	96.24
Deuterium Oxide	4.80 (DSS) 4.81 (TSP)		NA	NA	4.8	1.11	3.81	101.42	78.5	20.03
N-Dimethyl-formamide-d ₇	8.03 (1) 2.92 (5) 2.75 (5)	1.9 1.9	163.15 (3) 34.89 (7) 29.76 (7)	29.4 21.0 21.1	3.5	1.03	-61	153	36.7	80.14
Dimethyl Sulfoxide-d ₆	2.50 (5)	1.9	39.51 (7)	21.0	3.3*	1.19	18.55	189	46.7	84.17
1,4-Dioxane-d ₈	3.53 (m)		66.66 (5)	21.9	2.4	1.13	11.8	101.1	2.2	96.16
Ethanol-d ₆	5.19 (1) 3.56 (1) 1.11(m)		56.96 (5) 17.31 (7)	22 19	5.3	0.89	-114.1	78.5	24.5	52.11
Methanol-d ₄	4.87 (1) 3.31 (5)	1.7	49.15 (7)	21.4	4.9	0.89	-97.8	64.7	32.7	36.07
Methylene Chloride-d ₂	5.32 (3)	1.1	54.00 (5)	27.2	1.5	1.35	-95	39.75	8.9	86.95
Pyridine-d ₅	8.74 (1) 7.58 (1) 7.22 (1)		150.35 (3) 135.91 (3) 123.87 (3)	27.5 24.5 25	5	1.05	-41.6	115.2-115.3	12.4	84.13
1,2,2-Tetrachloroethane-d ₂	6.0		73.78 (3)			1.62	-44	146.5	8.20	169.86
Tetrahydrofuran-d ₈	3.58 (1) 1.73 (1)		67.57 (5) 25.37 (5)	22.2 20.2	2.4-2.5	0.99	-108.5	66	7.6	80.16
Toluene-d ₈	7.09 (m) 7.00 (1) 6.98 (5) 2.09 (5)	2.3	137.86 (1) 129.24 (3) 128.33 (3) 125.49 (3) 20.4 (7)	23 24 24 19	0.4	0.94	-95	110.6	2.4	100.19
Trifluoroacetic Acid-d	11.50 (1)		164.2 (4) 116.6 (4)		11.5	1.49	-15.4	72.4		115.03
Trifluoroethanol-d ₃	5.02 (1) 3.88 (4x3)	2(9)	126.3 (4) 61.5 (4x5)	22	5	1.41	-43.5	74.05		103.06

M.J. O'Neil, P.E. Heckelman, C.B.Koch, K.J. Roman, The Merck Index, an Encyclopedia of Chemicals, Drugs, and Biologicals – Fourteenth Edition, Merck Co., Inc. Whitehouse Station, NJ 2006.

The ¹H spectra of the residual protons and ¹³C
Approximate values only, may vary with pH, spectra were obtained on a Varian Gemini 200 spectrometer at 29.5°K. The NMR solvents used to acquire these spectra contain a maximum of 0.05% and 1.0% TMS (v/v) respectively. Since deuterium has a spin of 1, triplets arising from coupling to deuterium have the intensity ratio of 1:1:1. "m" denotes a broad peak with some fine structures. It should be noted that chemical shifts can be dependent on solvent, concentration and temperature.

concentration and temperature.

Melting and boiling points are those of the corresponding unlabeled compound (except for D_2O). These temperature limits can be used as a guide to determine the useful liquid range of the solvents. Information gathered from the Merck Index – 14th Edition.

★ HOD Peaks – NMR spectra of "neat" deuterated solvent always exhibit a peak due to H_2O in addition to the residual solvent peak. When the exchange rate between H₂O and HDO is slow on the NMR timescale the water peak appears as two peaks, a singlet corresponding to H₂O and a 1:1:1 triplet corresponding to HDO.